For Research Use Only. Not for use in diagnostic

c-Myc-tagged Protein PURIFICATION CARTRIDGE

(Clone 1G4)

CODE No. 3306K

*PURIFICATION to maintain protein activity
from eukaryote cell lysate and culture supernatant*

Product Description

Recombinant tagged proteins are widely used for protein purification and protein characterization. There are many purification methods and kits available. c-Myc epitope tag (EQKLISEEDL) has been widely used as a multi-purpose tag for protein purification.

MBL's c-Myc-tagged Protein PURIFICATION CARTRIDGE is ready-to-use-cartridge for the isolation of c-Myc-tagged protein from cell culture supernatants and cell lysate under neutral pH condition. The cartridge can be used directly on automated ÄKTA FPLC Systems.

The elution of c-Myc-tagged proteins from the cartridge is achieved by the addition of the c-Myc-tag peptide (EQKLISEEDL). As the c-Myc-tag peptide competes with c-Myc-tagged proteins on the cartridge, the purified c-Myc-tagged proteins maintain function and retain protein activity.

Components

	Quantity
	CODE No. 3306K
<u>c-Myc-tagged Protein PURIFICATION CARTRIDGE</u>	<u>1 mL × 1</u>
• c-Myc-tag cartridge (1/16 inch female end) × 1, Stop plug (1/16 inch male) × 2	

Elution Peptide (EQKLISEEDL) is NOT included.

It is available as CODE No. 3300-205 (1 mg × 5 vials).

Storage

Store for up to 1 year from date of receipt at 2-8°C. Do not freeze.

Cartridge Specifications

Specifications	
Column volume	1 mL
Cartridge body material	Polypropylene
Connector size	1/16 inch
Matrix	Cross-linked agarose
Ligand	Anti-c-Myc-tag mAb (clone: 1G4)
Binding capacity	Approx. 0.3 mg recombinant c-Myc-tagged Protein (32 kDa) /mL gel
Maximum pressure	0.3 MPa
Recommended flow rate	0.5-1 mL/min.
Storage buffer	0.1% ProClin 150/PBS

Material Preparation

Prepare the following reagents before affinity purification.

1. Lysis buffer : Suitable Lysis buffer varies with a kind of the c-Myc-tagged protein
(See **Additional Information**).
[Example of buffer constitution]
10-50 mM Tris-HCl (pH 7.5)
100-300 mM NaCl
1% NP-40 or Triton X-100
If necessary add Protease Inhibitor Cocktail
(e.g. SIGMA: code P8340, PIERCE: code 78415)
2. Washing buffer : Suitable Washing buffer varies with a kind of the c-Myc-tagged protein
(See **Additional Information**).
[Example of buffer constitution]
0.05% Tween-20 /PBS
To prevent non-specific protein binding to the gel,
0.05% Tween-20 should be added to Washing buffer.
3. Elution buffer : 0.5 mg/mL c-Myc-tag peptide in PBS.
4. Regeneration buffer : 0.17 M Glycine-HCl (pH 2.3)
5. Column storage buffer : 0.1% ProClin 150/PBS
6. A suitable liquid chromatography system with 1/16 inch tubing (e.g. GE Healthcare: ÄKTA system)

Note: For best results, filter all the buffers through 0.45 µm filters and degas before use.

Protocols

A. Sample preparation

The cartridge is optimized under only native conditions. It is not recommended for use under denaturing conditions or for purification of aggregated, unstable, and insoluble protein (e.g. inclusion bodies). Proteins solubilized with such as 6 M Guanidine-HCl or 8 M Urea cannot be purified using this cartridge (see **Additional Information**).

Cellular debris and particulate matter must be removed prior to purification. The protein extract should be centrifuged (10,000-20,000 × g for 15 min.) and filtered with a 0.45 µm filter to remove any remaining cells and particulates.

Highly viscous samples containing chromosomal DNA or RNA should be sonicated or treated with nuclease to reduce viscosity.

B. Purification

1. Fill the pump A and B in liquid chromatography system with Washing buffer and Elution buffer.
2. Remove one end of the stop plug from the cartridge, and start pumping Washing buffer at a flow rate of 0.5 mL/min. until a few drops fill in the top of the cartridge. Connect the cartridge to the pump side tubing.
3. Immediately remove the other end of the stop plug from the cartridge, connect the cartridge to the UV monitor side tubing.
4. Equilibrate the cartridge with 5-10 bed volumes of the Washing buffer at a flow rate of 1 mL/min.
5. Apply the sample to the cartridge at a flow rate of 0.5-1 mL/min. Collect fractions.
6. Wash the cartridge with 20 bed volumes of the Washing buffer until the absorbance approaches baseline at a flow rate of 1 mL/min.
7. Elute with 8 bed volumes of Elution buffer at a flow rate of 0.5-1 mL/min., and collect 0.5 mL fractions.

C. Regeneration and storage

1. Wash the cartridge with 10 bed volumes of Regeneration buffer.
2. Immediately wash the column with 10 bed volumes of Column storage buffer.
3. Screw the stop plug tightly, store the cartridge at 2-8°C.



Note: This cartridge may be used multiple times, depending on the usage conditions.

Related Products:

Please visit our website at <https://ruo.mbl.co.jp/>.

Example of Purification Result

Purification of N-terminus c-Myc-tagged GST

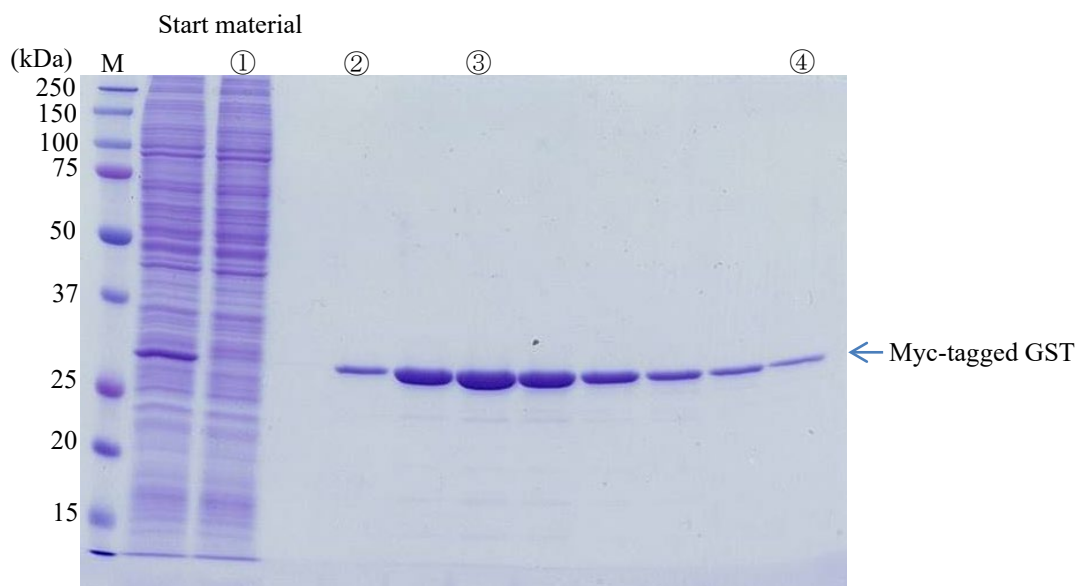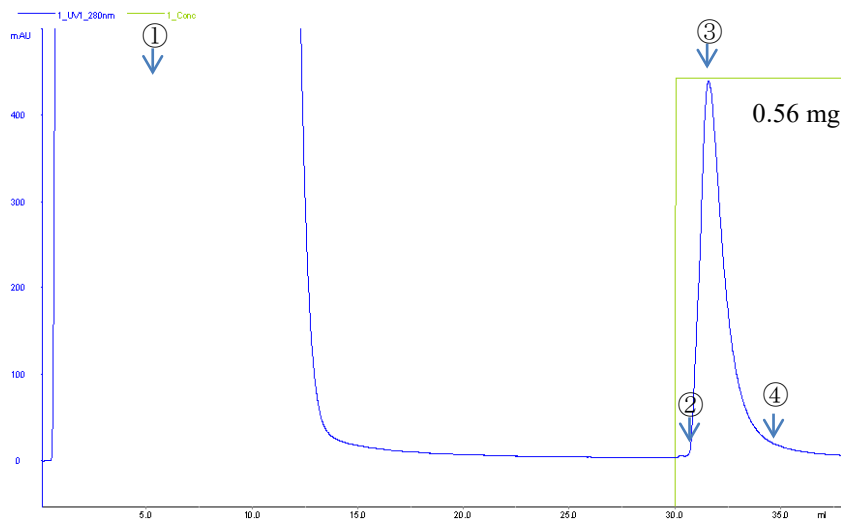Cartridge	: c-Myc-tagged Protein PURIFICATION CARTRIDGE
Sample	: c-Myc-tagged GST/293T (5×10^7 cells)
Cell lysis buffer	: 10 mM Tris-HCl, 150 mM NaCl, 1% NP-40 (pH 7.5)
Wash buffer	: 0.05% Tween-20/PBS (pH 7.2)
Elution buffer	: 0.5 mg/mL c-Myc peptide in PBS
Flow rate	: 0.75 mL/min.
Chromatography system	: ÄKTAexplorer 10S (GE Healthcare)

Additional Information

Several reagents were examined whether or not they were suitable for use with the c-Myc-tagged Protein PURIFICATION CARTRIDGE. The results are listed below.

Chaotropic agents

Urea	1 M	Yes
Guanidine-HCl	1 M	No

Reducing agents

DTT	10 mM	Yes
2-Mercaptoethanol	10 mM	Yes

Surfactants

Nonionic	Tween-20	5%	Yes
	TritonX-100	5%	Yes
	NP40	5%	Yes
	Digitonin	1%	Yes
	n-Octyl-β-D-gulcoside	1%	Yes
Zwitterionic	CHAPS	1%	Yes
	CHAPSO	1%	Yes
Anionic	SDS	0.05%	No
	Sodium Deoxycholate	0.1%	Yes
	Sodium Deoxycholate	0.5%	No

Others

NaCl	1 M	Yes
Glycerol	10%	Yes
EDTA	10 mM	Yes

The “Yes” indicates the reagents can be used in the Lysis buffer for this cartridge up to the indicated concentration. The “No” indicates the reagents cannot be used in the Lysis buffer for this cartridge at the indicated concentration.

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